

RapidScan® Cassette and Test Strip Orientation

ABRAXIS® Algal Toxin and Glyphosate test strips are compatible with and have been validated for use with the RapidScan ST-5 Reader. To run a test on the RapidScan® Test Strip Reader, the test strip and cassette must be oriented such that the camera within the reader is able to see both the test and control lines. This strip/cassette orientation permits it to determine the line intensities and provide a result.

The test strip should be placed in the cassette such that the wick end of the test strip is facing away from the numbered end of the cassette (arrows should be pointing away from the numbered end), as shown in the image below.



The cassette is then placed into the reader such that the numbered end of the cassette faces out of the reader, as shown in the images below.



Gold Standard Diagnostics Horsham, Inc.

124 Railroad Drive, Warminster, PA 18974 USA

(215) 357-3911 | info.abraxis@us.goldstandarddiagnostics.com | Website: www.abraxiskits.com

Then, interpret the test strip results following the instructions in the User's manual.

For ordering or technical assistance contact:

Gold Standard Diagnostics

124 Railroad Drive

Warminster, PA 18974

WEB: www.abraxiskits.com

Phone: (215) 357 3911

Fax: (215) 357 5232

Ordering: info.abraxis@us.goldstandarddiagnostics.com

Technical Support: support.abraxis@us.goldstandarddiagnostics.com

Date this Technical Bulletin is effective: 20SEP2023

Version: 02